

## Micro Mate

## l ow Volume Powder Sampler

The MicroMate is designed for low volume sampling of free flowing powders and granules.


The MicroMate can be easily adjusted so that it can take a range of sample sizes.

- Made from high quality 316 stainless steel
- Adjustable takes a range of sizes
- Simple to use



## Operation

- Insert the sampler into the product. Ensure that the inner rod extends beyond the outer tube.
- 2. At the required depth pull up the inner rod to draw the product into the sampler
- 3. Withdraw sampler.
- 4. Push down handle to release sample

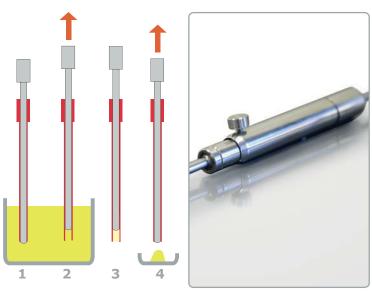





| Stock No.    | Sample Volume          | Length | Material | <b>Outer Diameter</b> |
|--------------|------------------------|--------|----------|-----------------------|
| 675 1070A350 | Variable - up to 0.1ml | 350mm  | 316ss    | 5mm                   |
| 675 1070A500 | Variable - up to 0.1ml | 500mm  | 316ss    | 5mm                   |

QUICK QUOTE ONLINE PRICING