



**QAQC LAB**

[www.qclabequipment.com](http://www.qclabequipment.com)

## Powder Thief

Best Selling Powder Thief


The PowderThief was originated over 20 years ago and is now a world leader in powder sampling. It is extensively used throughout the pharmaceutical, chemical and food industries around the world.

The PowderThief is suitable for sampling free flowing powders and granules. A range of tips is available so that a single sampler body can be used to accurately sample a range of different volumes.

- Made from high quality 316 stainless steel
- Wide range of sizes
- Simple to use




Tips - various sizes can be attached to the same body to allow different volumes to be sampled

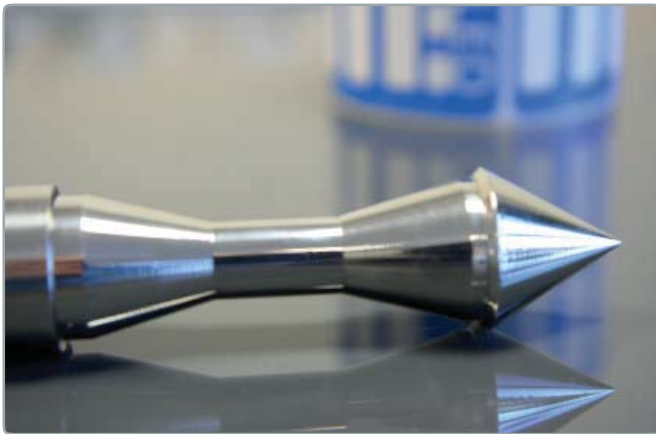

### Operation

1. Insert the sampler into the product, ensure that the tip is inside the sampler body.
2. At the required depth pull up the body to expose the tip. Powder will flow in around the tip.
3. Push down body of the sampler to trap the sample.
4. Withdraw sampler.
5. Pull up body to release the sample.

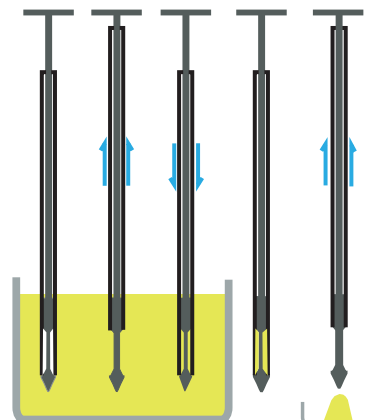



PowderThief samplers are available in a range of different sizes

### Sampling Tips

A range of standard tip sizes is available for each body diameter. Special sizes available on request.

Stock No.	Length	Diameter
675 1030A300	300mm	12.5mm
675 1030A600	600mm	12.5mm
675 1030A1000	1000mm	12.5mm
675 1030A1500	1500mm	12.5mm
675 1030A2000	2000mm	12.5mm
675 1030B600	600mm	19mm
675 1030B1000	1000mm	19mm
675 1030B1500	1500mm	19mm
675 1030B2000	2000mm	19mm
675 1030C600	600mm	25mm
675 1030C1000	1000mm	25mm
675 1030C1500	1500mm	25mm
675 1030D600	600mm	32mm
675 1030D1000	1000mm	32mm
675 1030D1500	1500mm	32mm
675 1030D2000	2000mm	32mm


Stock No.	Vol.	Diameter
675 1030AT025	0.25ml	12.5mm
675 1030AT050	0.50ml	12.5mm
675 1030AT075	0.75ml	12.5mm
675 1030AT100	1.00ml	12.5mm
675 1030BT2	2ml	19mm
675 1030BT3	3ml	19mm
675 1030BT5	5ml	19mm
675 1030BT8	8ml	19mm
675 1030BT10	10ml	19mm
675 1030CT10	10ml	25mm
675 1030CT20	20ml	25mm
675 1030CT25	25ml	25mm
675 1030DT20	20ml	32mm
675 1030DT25	25ml	32mm
675 1030DT50	50ml	32mm
675 1030DT100	100ml	32mm


[QUICK QUOTE](#)

[ONLINE PRICING](#)

... also available is a single use, disposable PowderThief


The Disposable PowderThief works in the same way as the stainless steel Powder Thief (above). Disposable samplers have the advantage that they can be used straight from the bag and never need to be cleaned, thereby saving you time and money.

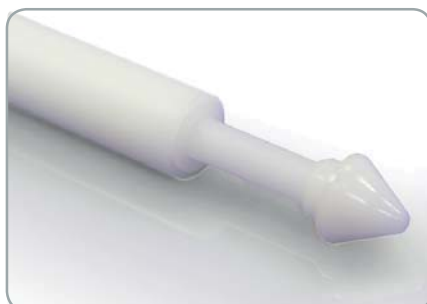

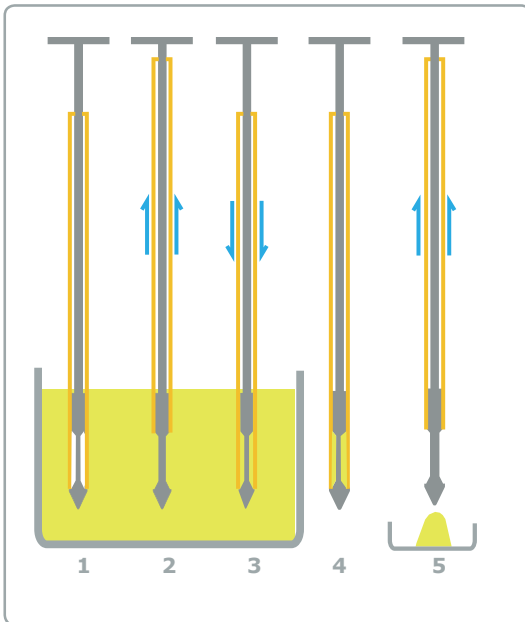

#### Operation

1. Insert the sampler into the product. Ensure that the tip is inside the sampler body
2. At the required depth pull up the body to expose the tip. The sample will fall in around the tip
3. Push down body of the sampler to trap the sample
4. Withdraw sampler
5. Pull up body to empty sample

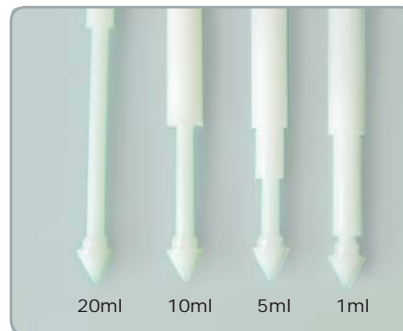

Available with a range of different tip sizes: 1ml, 5ml, 10ml and 20ml

Stock No.	Pre-Sterile	Capacity	Length	Diameter	Material	Per Box
675 8060H501	No	1ml	500mm	21mm	HDPE	20
675 8060H501S	Yes	1ml	500mm	21mm	HDPE	20
675 8060H505	No	5ml	500mm	21mm	HDPE	20
675 8060H505S	Yes	5ml	500mm	21mm	HDPE	20
675 8060H510	No	10ml	500mm	21mm	HDPE	20
675 8060H510S	Yes	10ml	500mm	21mm	HDPE	20
675 8060H520	No	20ml	500mm	21mm	HDPE	20
675 8060H520S	Yes	20ml	500mm	21mm	HDPE	20
675 8060H1001	No	1ml	1000mm	21mm	HDPE	20
675 8060H1001S	Yes	1ml	1000mm	21mm	HDPE	20
675 8060H1005	No	5ml	1000mm	21mm	HDPE	20
675 8060H1005S	Yes	5ml	1000mm	21mm	HDPE	20
675 8060H1010	No	10ml	1000mm	21mm	HDPE	20
675 8060H1010S	Yes	10ml	1000mm	21mm	HDPE	20
675 8060H1020	No	20ml	1000mm	21mm	HDPE	20
675 8060H1020S	Yes	20ml	1000mm	21mm	HDPE	20

**QUICK QUOTE**

**ONLINE PRICING**

Note: Other tip sizes are available between 1 and 20ml - subject to minimum order quantities  
1500mm long samplers are available on request - subject to minimum quantities